



# Chemical Compatibility of UF Products

## Chemical Compatibility of UF Products

### Solvents, Acids, Bases

Key to the above table:

R = Recommended

NR = Not Recommended

TST = Testing prior to use is recommended

ND = No Data currently available. Under investigation

#### SOLVENTS

	Acetone	Acetonitrile	Acetonitrile (10%)	Acetonitrile (40% in 1% TFA)	Alconox (1%)	Ammonium Acetate	Ammonium Sulfate (50%)	Ammonium Sulfate	Amyl Alcohol	Benzyl Alcohol (1%)	n-Butanol	CAPS (250mM, pH11)	CHAPS (100mM)	Deco 90 (1%)
Microcon	NR	NR	R	TST	R	R	R	R	NR	R	NR	R	R	R
Centricon	NR	NR	R	TST	R	R	R	R	NR	R	NR	R	R	R
Centriprep	NR	NR	R	TST	R	R	R	R	NR	R	NR	R	R	R
Centriplus	NR	R	R	R	R	R	R	R	NR	R	NR	R	R	R
Centricon Plus with Biomax-PB Membrane	NR	R	R	R	R	R	R	R	R	R	TST	R	R	R
Centricon Plus with Ultracel-PL Membrane	R	R	R	R	R	R	R	R	R	R	TST	R	R	R
Amicon Ultra	NR	ND	R	ND	ND	ND	R	R	ND	ND	ND	ND	ND	ND
Ultrafree MC/CL with PVDF Membrane	NR	R	R	TST	R	R	R	R	R	R	TST	TST	R	R
Ultrafree MC/CL with Biomax-PB Membrane	NR	R	R	TST	R	R	R	R	R	R	TST	R	R	R
Ultrafree MC/CL with Ultracel-PL Membrane	R	R	R	TST	R	R	R	R	R	R	TST	R	R	R
Ultracel YM/PL Membrane Disc	R	R	R	R	R	R	R	R	R	R	TST	R	R	R
Biomax-PB Membrane Disc	NR	R	R	R	R	R	R	R	R	R	TST	R	R	R

8000 Series Stirred Cells	NR	R	R	R	R	R	R	R	R	R	TST	R	R	R
2000 Series Stirred Cells	NR	NR	TST	NR	R	R	R	R	NR	R	NR	R	R	R

---

© 1994-2008 Millipore Corporation. All rights reserved.