




Data Sheet

VOC-Pak™ Point-of-Use Polisher

Final purification step at the ultrapure water point-of-delivery efficiently removes Volatile Organic Compounds (VOCs)

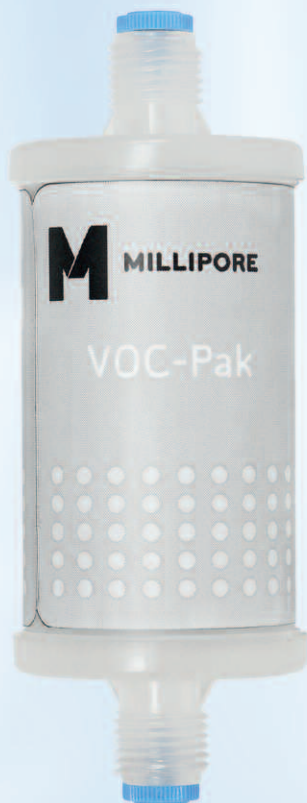

VOCs are industrial contaminants that may cause health safety issues and therefore should be detected in the environment. VOCs are routinely analyzed in water, soil, soil gas and air.

Specifications for VOCs in drinking water have been established by different regulatory agencies in many countries. A table showing the Maximum Allowable Concentration (MAC) for some of the different VOCs monitored in the United States, Europe, Japan and China, as well as by the World Health Organization (WHO), is available on the Millipore web site:


www.millipore.com/labwater

Benefits

- The VOC-Pak polisher is easily connected to the outlet of all Millipore Type I water purification systems (Milli-Q®, Direct-Q®, Synergy® and Simplicity® systems)
- The final purification step provides VOC-free ultrapure water at high flow rate—when you need it
- Delivers a minimum of 300 l of VOC-free ultrapure water
- Validated for efficient removal of VOCs
- Each VOC-Pak cartridge is delivered with a Certificate of Quality

VOC-PAK POLISHER DESIGN

The VOC-Pak cartridge is designed to assure scientists performing VOC analysis that the ultrapure water they use to rinse glassware, produce standard solutions or dilute samples is VOC-free.

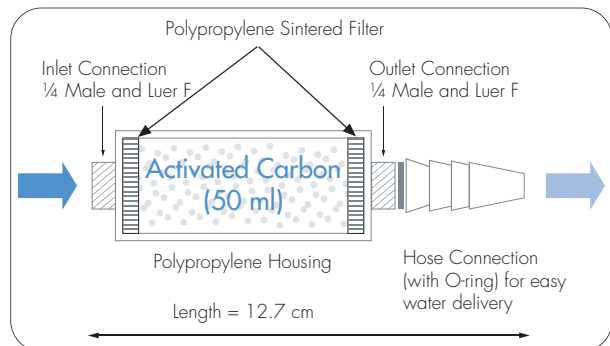

The purification media used to achieve this goal is a specific activated carbon selected after tests performed on media from many suppliers.

VOC-PAK POLISHER OPERATION

The VOC-Pak cartridge is easy to use: simply connect the VOC-Pak unit to the outlet of the

Millipore ultrapure water system and flush 5 l of water. You are now ready to deliver at least 300 l of VOC-free ultrapure water.

VOC-PAK POLISHER SPECIFICATIONS

The VOC-Pak cartridge is designed to produce water in which the VOC concentration is below the analytical quantification limit (**) for the VOCs listed in Table 1, provided that the VOC concentration in feed water does not exceed 1 ppb (1 µg/l), and the Total Oxidizable Carbon (TOC) of the ultrapure water entering the VOC-Pak cartridge is below 5 ppb.

In order to warrant this water quality, the VOC-Pak cartridge was qualified by a challenge test: the VOC-Pak cartridge was fed by ultrapure water containing a 1 ppb (1 µg/l) concentration of the VOC to be eliminated.

The concentration of the VOC was measured by Gas Chromatography (GC) downstream of the VOC-Pak cartridge after passage of 50, 100, 150, 200, 250, 300 and 350 l in order to determine the breakthrough point of the different VOCs.

The results are displayed in Table 1 for a processed volume of 350 l.

Table 1:

Results of the VOC-Pak challenge test for a processed volume of 350 l

| VOC tested | Challenge (ppb) (*) | Quantification limit (ppb) (**) | Concentration after VOC-Pak (ppb) (***) |
|---------------------------|---------------------|---------------------------------|---|
| benzene | 0.97 | < 0.05 | < 0.05 |
| bromodichloromethane | 1.56 | < 0.05 | < 0.05 |
| bromoform | 1.58 | < 0.50 | < 0.50 |
| carbon tetrachloride | 1.06 | < 0.05 | < 0.05 |
| chlorobenzene | 0.94 | < 0.05 | < 0.05 |
| chloroform | 1.07 | < 0.05 | < 0.05 |
| dibromochloromethane | 1.01 | < 0.10 | < 0.10 |
| 1,2-dichlorobenzene | 1.01 | < 0.05 | < 0.05 |
| 1,4-dichlorobenzene | 1.53 | < 0.05 | < 0.05 |
| 1,2-dichloroethane | 1.08 | < 0.05 | < 0.05 |
| 1,1-dichloroethene | 0.87 | < 0.05 | < 0.05 |
| cis-1,2-dichloroethene | 0.95 | < 0.05 | < 0.05 |
| trans-1,2-dichloroethene | 1.53 | < 0.05 | < 0.05 |
| 1,2-dichloropropane | 0.86 | < 0.05 | < 0.05 |
| cis-1,3-dichloropropene | 0.97 | < 0.04 | < 0.04 |
| trans-1,3-dichloropropene | 0.91 | < 0.05 | < 0.05 |
| ethylbenzene | 1.09 | < 0.05 | < 0.05 |
| tetrachloroethene | 1.44 | < 0.05 | < 0.05 |
| toluene | 0.96 | < 0.10 | < 0.10 |
| 1,1,1-trichloroethane | 0.94 | < 0.10 | < 0.10 |
| 1,1,2-trichloroethane | 0.93 | < 0.05 | < 0.05 |
| trichloroethene | 1.02 | < 0.05 | < 0.05 |
| o-xylene | 0.79 | < 0.05 | < 0.05 |
| m-xylene | 1.02 | < 0.05 | < 0.05 |
| p-xylene | 0.94 | < 0.05 | < 0.05 |

(*) Challenge = VOC concentration in feed water upstream of VOC-Pak cartridge

(**) Quantification limit: limit of quantification of the VOC measured by GC

(***) Concentration of VOC below the VOC-Pak cartridge for a feed water concentration of 1 ppb



Q-POD dispenser fitted with a VOC-Pak cartridge

Analysis of the VOCs in ultrapure water was performed according to Millipore procedure 8421450, using an Eclipse 4660 Purge-and-Trap sample concentrator from OI Corporation™ (College Station, Texas) and a GC-MS system (GC model 6890N and MS model 5975B) from Agilent Technologies®, Inc., Santa Clara, California. A copy of the complete qualification report with

description of the analytical procedure is available upon request.


Additionally, each lot of VOC-Pak cartridges is analyzed to ensure that the following VOCs (Table 2) are below the analytical quantification limit in water produced by a VOC-Pak fed with ultrapure water from a Millipore Milli-Q system (resistivity at 18.2 MΩ·cm at 25 °C and TOC < 5 ppb).

Table 2:

Quantification limits (ppb) of VOCs analyzed in each lot of VOC-Pak cartridges

| Compound | QL* (ppb) | Compound | QL (ppb) | Compound | QL (ppb) |
|------------------------------------|-----------|--------------------------------------|----------|----------------------------|----------|
| benzene | < 0.05 | 1,3-dichlorobenzene | < 0.05 | styrene | < 0.10 |
| bromobenzene | < 0.10 | 1,4-dichlorobenzene | < 0.05 | 1,1,1,2-tetrachloroethane | < 0.05 |
| bromochloromethane | < 0.05 | 1,1-dichloroethane | < 0.05 | 1,1,2,2-tetrachloroethane | < 0.10 |
| bromodichloromethane | < 0.05 | 1,2-dichloroethane | < 0.05 | tetrachloroethene | < 0.05 |
| bromoform | < 0.50 | 1,1-dichloroethene | < 0.05 | toluene | < 0.10 |
| bromomethane | < 0.20 | cis-1,2-dichloroethene | < 0.05 | 1,2,3-trichlorobenzene | < 0.05 |
| n-butylbenzene | < 0.05 | trans-1,2-dichloroethene | < 0.05 | 1,2,4-trichlorobenzene | < 0.05 |
| sec-butylbenzene | < 0.50 | 1,2-dichloropropane | < 0.05 | 1,1,1-trichloroethane | < 0.10 |
| tert-butylbenzene | < 0.05 | 1,3-dichloropropane | < 0.10 | 1,1,2-trichloroethane | < 0.05 |
| carbon tetrachloride | < 0.05 | 2,2-dichloropropane | < 0.05 | trichloroethene | < 0.05 |
| chlorobenzene | < 0.05 | 1,1-dichloropropene | < 0.05 | trichloromonofluoromethane | < 0.10 |
| chloroethane | < 1.00 | cis-1,3-dichloropropene | < 0.05 | 1,2,3-trichloropropane | < 0.05 |
| chloroform | < 0.05 | trans-1,3-dichloropropene | < 0.05 | 1,2,4-trimethylbenzene | < 0.05 |
| 2-chlorotoluene | < 0.05 | ethylbenzene | < 0.05 | 1,3,5-trimethylbenzene | < 0.10 |
| 4-chlorotoluene | < 0.10 | hexachlorobutadiene | < 0.05 | o-xylene | < 0.05 |
| dibromochloromethane | < 0.10 | isopropylbenzene | < 0.05 | m-xylene | < 0.05 |
| 1,2-dibromo-3-chloropropane (DBCP) | < 0.05 | 4-isopropyltoluene | < 0.05 | p-xylene | < 0.05 |
| 1,2-dibromoethane | < 0.05 | methylene chloride (dichloromethane) | < 0.10 | vinyl chloride | < 0.20 |
| dibromomethane | < 0.10 | naphtalene | < 0.10 | | |
| 1,2-dichlorobenzene | < 0.05 | n-propylbenzene | < 0.50 | | |

*QL = Quantification Limit – expressed in ppb (µg/l)

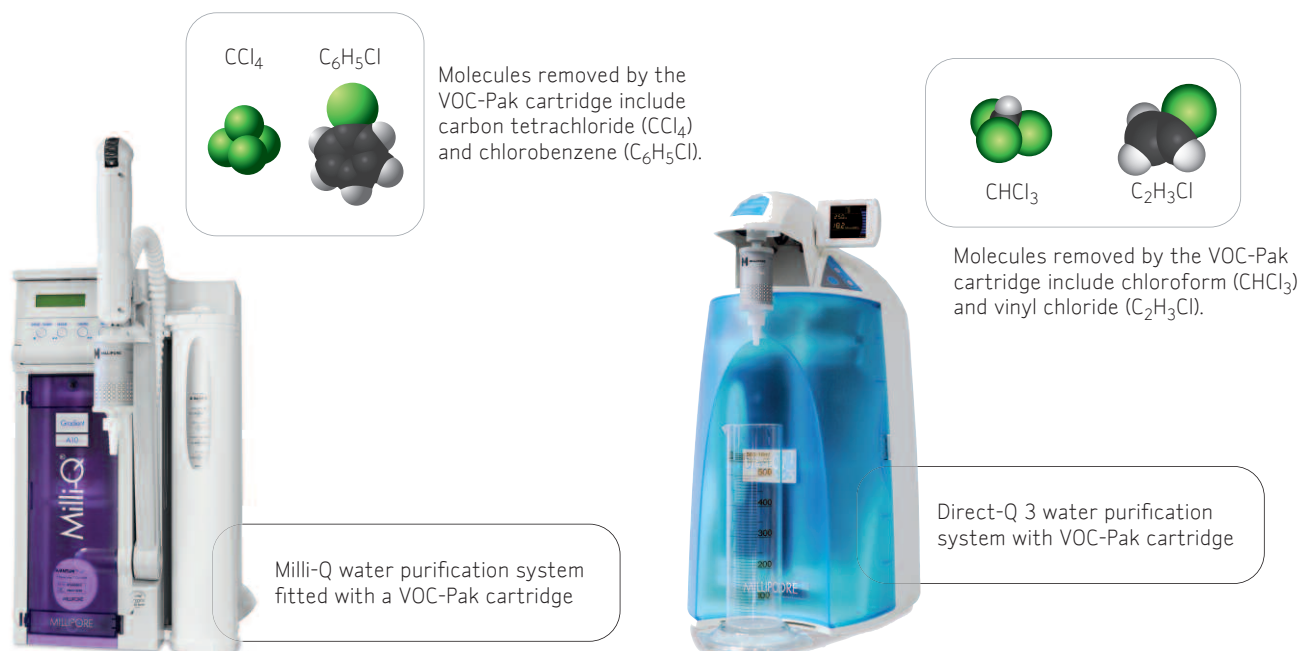

ORDERING INFORMATION

| Description | Catalogue Number |
|--|------------------|
| VOC-Pak cartridge (1/pk) delivered hermetically sealed, with a Certificate of Quality | VOCPAK001 |
| Cartridge installation and Conditioning Kit including the following reusable parts: | EDSKIT001 |
| <ul style="list-style-type: none">• Polyethylene 1/4 Gaz F-Hose barb connection with O-ring• Polyethylene 1/4 Gaz F – 1/4 Gaz F connectors for connection to Millipak® final filter | |

Millipore offers more innovative technologies and stronger application support to streamline processes and provide consistently reliable results. Our Lab Water experts take the time to evaluate the needs of individual labs and particular applications in order to recommend a system that balances water quality with volume and distribution requirements, removing water quality concerns so customers can focus on their research.



www.millipore.com/offices

ADVANCING LIFE SCIENCE TOGETHER™
Research. Development. Production.

Millipore, Milli-Q, Direct-Q, Synergy, Simplicity, and Millipak are registered trademarks of Millipore Corporation.
VOC-Pak, the M mark and Advancing Life Science Together are trademarks of Millipore Corporation.
OI Corporation is a trademark of OI Corporation. Agilent Technologies is a registered trademark of Agilent Technologies, Inc.
Lit. No. PB1632EN00 Printed in France 04/09
© 2009 Millipore Corporation, Billerica, MA 01821 U.S.A.
All rights reserved. Design: Un Coin de Paradis