

Aerex™ 2 Hydrophobic Cartridge Filters

■ High performance gas filters for use in industrial fermentation applications



Aerex 2 0.2 µm hydrophobic cartridge filters provide high flow rates at low pressure differentials. Designed to remove bacteria and viruses from moist gas streams, the filters are ideally suited for use in industrial fermentation applications. Aerex 2 cartridge filters are manufactured using PTFE membrane and are 100 % integrity tested to assure optimal sterilizing performance. Each filter lot also must pass rigorous stress testing before release.

Cost Effective

High performance Aerex 2 cartridge filters offer significant savings of operating costs and capital. The combination of high flow rates and low pressure differentials lowers the air compressor loads that are required to achieve the desired flows. This reduction lowers the energy costs of air compressor operation as well as the required frequency of air compressor maintenance. The efficiency of Aerex 2 filters also allows the use of smaller housings, thereby reducing the required capital costs.

Advantages Provided by Aerex 2

- Sterility Assurance
- Improved Process Efficiency
- Long Service Life
- Quality Assurance

Sterility Assurance

Proven Bacterial and Viral Retention

Millipore's qualification of the AereX 2 cartridge filter includes a challenge regimen designed to test viral and bacterial retention by two challenge mechanisms: aerosol and grow-through. Challenge tests are completed on filters that have been exposed to 200 steam-in-place cycles. This challenge test protocol demonstrates organism retention over a long service life, not just when the filter is new.

Bacterial Growththrough Resistance

Millipore's retention test simulates the service challenge of retaining bacteria even after long use. The test entails first exposing AereX 2 filter cartridges to 200 thirty-minute steam-sterilization cycles in the forward direction at 145 °C. The test continues with the exposure of the cartridges to a moist gas stream containing aerosolized bacteria in a concentration $>1 \times 10^7$ cfu/cm². Under these rigorous test conditions, AereX 2 retained all *Brevundimonas diminuta* for 21 days.

Virus (Bacteriophage) Retention

Millipore's bacteriophage retention test performed after steam sterilization simulates the service challenge of retaining virus after long use. The test entails first exposing AereX 2 cartridge filters to 200 thirty-minute steam sterilization cycles in the forward direction at 145 °C. The test continues with the exposure of the filters to a moist gas stream containing the bacteriophage Φ X-174 at a concentration of 10^7 — 10^{10} virus particles per 10-inch cartridge. (The model bacteriophage Φ X-174 is icosahedral, has no tail required for infectivity, and is 29 nm in diameter.) These conditions represent a worst-case service challenge. After these exposures, the cartridges were tested for viral retention. AereX 2 met this challenge, retaining all bacteriophage particles under these extreme conditions.

Improved Process Efficiency

High Flow Rates

AereX 2 filters redesigned core allows for an open structure without compromising strength. This open structure provides superior flow rates, which minimizes the number of filter elements required to achieve a given flow volume, thereby increasing process efficiency. This process efficiency permits the use of smaller cartridge housings and reduces capital costs.

Low Pressure Differentials

AereX 2 filters operates at full efficiency without requiring the high pressure differentials usually needed to achieve high flow volumes. Low pressure differentials reduce air compressor loads, thereby saving energy. Low pressure differentials also preserve filters and air compressors, thereby reducing the need for compressor maintenance.

Long Service Life

The physical stress resulting from high pressure differentials and exposure to steam sterilization can result in the premature failure of hydrophobic

cartridge filters. AereX 2 filters are designed to withstand each of these challenges.

Validated Steam Sterilization


AereX 2 filters have been validated to withstand 200 steam-in-place cycles at 145°C. AereX 2 filters have maintained their integrity even after exposure to 200 steam-in-place cycles.


Validated Thermal and Hydraulic Stress Resistance

Millipore has demonstrated the ability of AereX 2 cartridges to resist hydraulically induced failure by validating their integrity against pulsed hydraulic pressure at 4.1 barg (60 psid) in the forward direction and 1.4 barg (20 psid) in the reverse direction.

Withstands Thermal Stress

AereX 2 filters have proved resistant to high levels of thermal stress. The multiple steam-in-place test is designed to demonstrate product resistance to thermally-induced stress. The test entails steaming the cartridges in place for 200 cycles at 145 °C. The test continues as cartridges are integrity tested after 200 steam cycles.

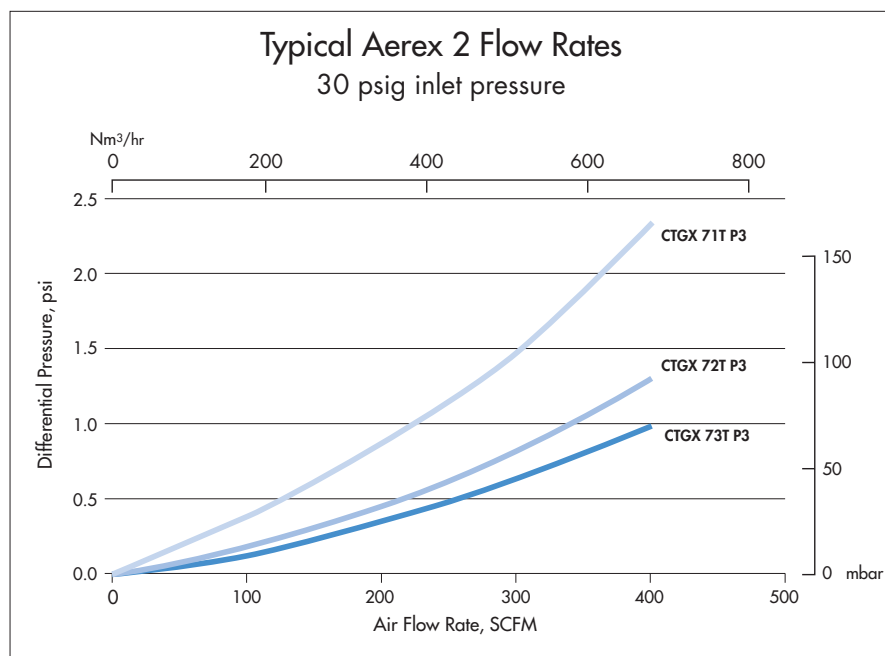

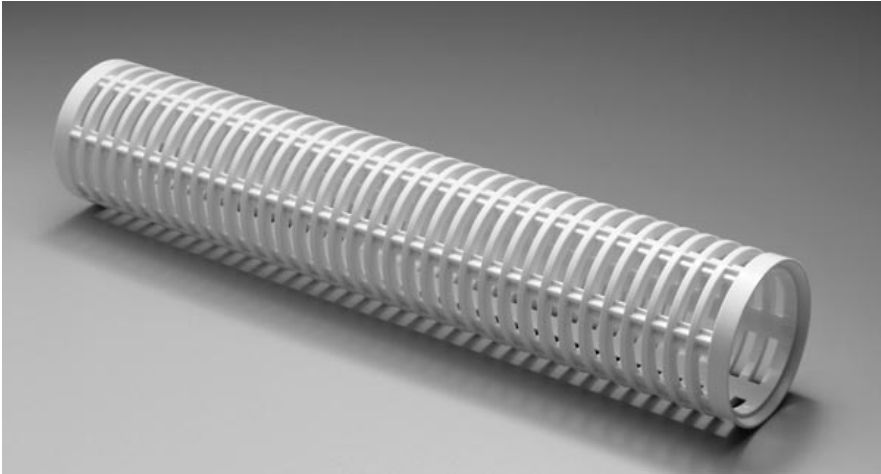

Aerex 2 filters open structured core provides superior flow rates.

Quality Assurance

A Certificate of Quality is provided with every Aerex 2 filter. These certificates record all relevant manufacturing lot-release criteria and accurately record quality test results, ensuring that each filter meets the minimum performance claims as detailed in the certificate. Millipore's Aerex 2 cartridges meet or surpass these claims. Aerex 2 cartridges are designed, developed, and manufactured in accordance with a Quality Management System that meets an ISO® 9000 Quality Systems Standard, as determined by an accredited registering body.

Specifications

Materials	PTFE membrane Spun-bonded polypropylene pleat supports upstream and downstream Polysulfone core Rigid polypropylene outer sleeve, end caps and adapter Silicone O-rings
Connections	Double 2-226 O-rings, bayonet with spear
Pore Size	0.2 µm
Nominal Cartridge Length	10 inch (25 cm) 20 inch (50 cm) 30 inch (75 cm)
Maximum Differential Pressure	<p><i>Forward Pressure:</i> 4.1 bar (60 psid) at 25 °C 340 mbar (5 psid) at 145 °C</p> <p><i>Reverse Pressure:</i> 1.4 bar (20 psid) at 25 °C (Intermittent)</p> <p>Aerex 2 hydrophobic cartridge filters must pass a hydraulic stress test prior to lot release.</p>
Effective Filtration Area	Approximately 0.66 m ² (7.1 ft ²) per 10-inch (25 cm) cartridge
Bacterial Growththrough Resistance	Aerex 2 cartridges resist the grow-through of aerosolized <i>B. diminuta</i> at a concentration of > 1x10 ⁷ cfu/cm ² for 21 days after 200 thirty-minute steam sterilization cycles at 145 °C.
Virus (Bacteriophage) Aerosol Retention	Retention of the model bacteriophage Φ X-174 was complete at exposure levels of 10 ⁷ — 10 ¹⁰ virus particles at 50 scfm (85 Nm ³ /hr) after 200 thirty-minute steam sterilization cycles at 145 °C. This rate of retention under these conditions represents a worst-case challenge for hydrophobic filters meeting these specifications.



Ordering Information

Description		Catalogue No.
Aerex 2, 10" Cartridge	3/pkg	CTGX 71T P3
Aerex 2, 20" Cartridge	3/pkg	CTGX 72T P3
Aerex 2, 30" Cartridge	3/pkg	CTGX 73T P3

To Place an Order or Receive Technical Assistance

For additional information call your nearest Millipore office:

In the U.S. and Canada, call toll-free **1-800-MILLIPORE (1-800-645-5476)**

In the U.S., Canada and Puerto Rico, fax orders to **1-800-MILLIFX (1-800-645-5439)**

On the Internet

<http://www.millipore.com>

E-mail: tech_service@millipore.com

Millipore Worldwide

AUSTRALIA

Tel. 1 800 222 111
or (02) 9888 8999
Fax (02) 9878 0788

AUSTRIA

Tel. (01) 877-8926
Fax (01) 877-1654

BALTIC COUNTRIES

Tel. +358 9 804 5110
Fax +358 9 256 5660

BELGIUM AND LUXEMBOURG

Tel. +32 2 726 88 40
Fax +32 2 726 98 84

BRAZIL

Tel. (011) 5548-7011
Fax (011) 5548-7923

CANADA

Tel. 1-800-645-5476
Fax 1-800-645-5439

CHINA, PEOPLE'S REPUBLIC OF

Beijing:
Tel. (8610) 8519 1250
(8610) 6518 1058
Fax (8610) 8519 1255
Guangzhou:
Tel. (8620) 8755 4049
Fax (8620) 8752 0172
Hong Kong:
Tel. (852) 2803 9111
Fax (852) 2513 0313
Shanghai:
Tel. (8621) 5306 9100
Fax (8621) 5306 0838

CZECH REPUBLIC

Tel. 02-2051 3841
Fax 02-2051 4298

DENMARK

Tel. 70 10 00 23
Fax 70 10 13 14

EASTERN EUROPE, C.I.S., AFRICA, MIDDLE EAST AND GULF

Life Sciences Division:
Tel. +33 3 88 38 9536
Fax +33 3 88 38 9539
BioPharmaceutical Division:
Tel. +43 1 877-8926
Fax +43 1 877-1654
Laboratory Water Division:
Tel. +33 1 30 12 7000
Fax +33 1 30 12 7180

FINLAND

Tel. (09) 804 5110
Fax (09) 256 5660

FRANCE

Tel. 01 30 12 7000
Fax 01 30 12 7180

GERMANY

Tel. (06196) 494-0
Fax (06196) 43901

HUNGARY

Tel. 1-205 9784
Fax 1-205 9792

INDIA

Tel. (91) 80-839 46 57
Fax (91) 80-839 63 45

IRELAND

Life Sciences and Laboratory Water Divisions:
Tel. +44 1923 816375
Fax +44 1923 818297
BioPharmaceutical Division:
Tel. (021) 883 666
Fax (021) 883 048

ITALY

Vimodrone (Milano):
Tel. (02) 25.07.81
Fax (02) 26.50.324
Roma:
Tel. (06) 52.03.600
Fax (06) 52.95.735

JAPAN

Tel. (03) 5442-9711
Fax (03) 5442-9736 Life Sciences
9737 BioPharm.
9734 Lab Water

KOREA

Tel. (822) 551-0990
Fax (822) 551-0228

MALAYSIA

Tel. 603-757 1322
Fax 603-757 1711

MEXICO

Tel. (525) 576-9688
Fax (525) 576-8706

THE NETHERLANDS

Tel. 076-5022000
Fax 076-5022436

NORWAY

Tel. 22 67 82 53
Fax 22 66 04 60

POLAND

Tel. 22-669 12 25
22-663 70 31
Fax 22-663 70 33

PUERTO RICO

Tel. (787) 273-8495
Fax (787) 747-6553

SINGAPORE

Tel. (65) 842 1822
Fax (65) 842 4988

SPAIN AND PORTUGAL

Tel. +34 917 283 960
Fax +34 917 292 909

SWEDEN

Tel. 08-628 6960
Fax 08-628 6457

SWITZERLAND

Volketswil:
Tel. (01) 908-30-60
Fax (01) 908-30-80
Lausanne:
Tel. (021) 641 2550
Fax (021) 641 2551

TAIWAN

Taipei:
Tel. (886-2) 700-1742
Fax (886-2) 755-3267
Hsin Chu City:
Tel. (886-3) 571-0178
Fax (886-3) 572-9520

U.K.

Tel. 01923 816375
Fax 01923 818297

U.S.A.

Tel. (781) 533-6000
Fax (781) 533-3110

IN ALL OTHER COUNTRIES

Millipore Intertech (U.S.A.)
Tel. +1 (781) 533-8622
Fax +1 (781) 533-8630

MILLIPORE

Millipore is a registered trademark of Millipore Corporation.

Aerex is a trademark of Millipore Corporation.

ISO is a registered trademark of The International Organization for Standardization.

Lit. No. PF1080EN00 Rev. 7/01 01-159

Printed in U.S.A.

Copyright 2000, Millipore Corporation. All rights reserved.