



## **Re-blot Plus**

**Strong Antibody Stripping Solution**

**2504**

**FOR RESEARCH USE ONLY  
Not for use in diagnostic procedures.**

This page left blank intentionally.

---

## Application

The MILLIPORE Re-Blot Plus Western Blot Strong Antibody Stripping Solution is effective for removal of antibodies from Western blots that have been developed with chemiluminescence or radioactive iodine or other isotopes. It is not recommended for stripping colorimetric substrates (TMB, DAB, 4-chloronaphthol, etc.), as it is not possible to effectively remove substrates that precipitate at the reaction site.

The Re-Blot Plus Western Blot Strong Antibody Stripping Solution should be used only for qualitative purposes until it has been established by comparative blot analysis that stripping does not quantitatively affect a given antigen.

*For Research Use Only; Not for use in diagnostic procedures*

---

## Test Principle

Western blotting is a commonly used technique for studying protein function and localization. Typically, protein samples are electro-phoresed on SDS-PAGE and transferred to a membrane such as nitrocellulose or nylon, where they are probed with specific antibodies. Unlike nucleic acid based technologies, which allow reuse of Southern and Northern blots, it has been difficult to reuse Western blots.

Stripping and re-probing of Western blots offers several advantages:

- 1) Conservation of samples that are expensive or available only in limited quantities,
- 2) Analysis of a given blot using several different antibodies, e.g. subtype- or isoform-specific antibodies,
- 3) Re-analysis of anomalous results and confirmation with the same or a different antibody,
- 4) Correcting errors in incubation with the wrong antibody,
- 5) Cost savings in reagents and time by reusing the same blot.

While antigen and antibody-based immunoaffinity matrices, such as Sepharose conjugates, have been reused many times without compromising antigen-antibody reactivity, the need for pH extremes and chaotropic agents has precluded the application of these methods to Western blotting.

The MILLIPORE Re-Blot Plus Western Blot Strong Antibody Stripping Solution contains specially formulated solutions that quickly and effectively remove antibodies from Western blots without significantly affecting the immobilized proteins.

Advantages of the Re-Blot Plus Western Blot Strong Antibody Stripping Solution include:

- No pungent-smelling  $\beta$ -mercaptoethanol is contained in the Antibody Stripping Solution.
- Antibody stripping is done at room temperature. No heating of blots is required.
- Blots can be stripped of antibodies in approximately 15 minutes at room temperature.
- Blots may be reused in 25 minutes.

---

## Kit Components

1. Strong Antibody Stripping Solution (10x) - 2 containers / 25 mL each

---

## Materials Not Supplied

- Standard Blot or blot strips, on nitrocellulose or PVDF/nylon membrane.
- Blocking Solutions.
- Plastic Wrap, such as Saran Wrap, for storage of blots that will not be re-probed immediately.
- Distilled Water, for reagent dilution.
- Plastic Trays for incubation of blots or blot strips in stripping, washing and blocking solutions.
- Positive and Negative Stripping Controls. It is recommended that one new strip (not subjected to stripping solution) be included for comparison purposes.

---

## Storage

The Strong Antibody Stripping Solution should be stored at 2-8°C upon arrival. Product is stable for 3 to 6 months after receipt. If Antibody Stripping Solution crystallizes upon storage, it may be re-dissolved with gentle warming at 37°C before use.

**Note:** To prevent reagent degradation secure the cap tightly upon storage. Avoid extended exposure to air.

---

## Preparation of Reagents

### 1. Dilution of Antibody Stripping Solution

#### IMMEDIATELY BEFORE USE:

Dilute Strong Antibody Stripping Solution 10x with distilled water to obtain a 1x solution. If Antibody Stripping Solution contains crystals, warm gently at 37°C until crystals have dissolved completely. Prepare enough solution to allow free movement of strips or blots during incubation, typically 4 mL per strip or 20 mL per standard blot. Use the following chart for suggested volumes of stripping or blocking solution.

| # of Strips or Blots  | Amount of 10x Antibody Stripping Solution | Amount of Distilled Water | Resulting Amount of 1x Working Stripping Solution |
|-----------------------|---|---------------------------|---|
| 1 Strip               | 400 µL                                    | 3.6 mL                    | 4 mL  |
| 5 Strips              | 1 mL                                      | 9.0 mL                    | 10 mL   |
| 1 Blot<br>(7 x 10 cm) | 2 mL                                      | 18.0 mL                   | 20 mL   |

#### Selection of Reagent

*Re-Blot Plus* Strong Antibody Stripping Solution gives good results on both nitrocellulose and PVDF membranes. However, the *Re-Blot Plus* Mild Antibody Stripping Solution (MILLIPORE Catalog Number 2502) may perform better with certain proteins and antibodies.

---

## Assay Instructions

**Note:** *The blots or individual strips that are to be re-used should be prepared for stripping immediately after their first usage. If stripping cannot be performed right away, membranes can be wrapped in plastic wrap and stored moist in PBS at 4°C. DO NOT STORE BLOTS IN DRY FORM.*

1. Blot Stripping. Fill plastic tray with appropriate amount of 1x Antibody Stripping Solution (see Preparation of Reagents for suggested quantities).
2. Using tweezers or forceps, submerge blot or blot strips in stripping solution. Incubate with gentle mixing for 15 minutes at room temperature.

Note: It may be necessary to increase the stripping incubation time when using blots that have been stripped previously. Simply increase the stripping time by 5 to 10 minutes, if needed.

3. Blocking. Fill a clean plastic tray with an equal amount of blocking buffer. Conventional blocking buffers such as 20 mM Tris HCl, pH 8.0; 150 mM NaCl; 0.1% Tween 20; 5% dry milk or similar are appropriate.
4. Wash blots two times 5 minutes each with blocking buffer.
5. The blot is now ready for reprobing with antibodies. It is suggested that users employ their own proven protocol and chemiluminescent detection method.

---

## Technical Hints

To increase efficiency of the kit it is recommended to use minimal sufficient amounts of antigens transferred to the membranes. When several antigens are to be detected sequentially, it is recommended to start with antigens from which weaker signal is expected.

The proper use of the Re-Blot Plus Western Blot Strong Antibody Stripping Solution should not drastically affect membrane bound antigens. However, this kit should only be used for qualitative purposes, unless it has been established that stripping does not quantitatively affect a given antigen. Most antigens should withstand at least 5 stripping cycles. However, it is to be taken into consideration that during each stripping cycle small portions of membrane-immobilized proteins will be removed.

---

## Warranty

**Millipore Corporation** ("Millipore") warrants its products will meet their applicable published specifications when used in accordance with their applicable instructions for a period of one year from shipment of the products. **MILLIPORE MAKES NO OTHER WARRANTY, EXPRESSED OR IMPLIED. THERE IS NO WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE.** The warranty provided herein and the data, specifications and descriptions of Millipore products appearing in Millipore's published catalogues and product literature may not be altered except by express written agreement signed by an officer of Millipore. Representations, oral or written, which are inconsistent with this warranty or such publications are not authorized and if given, should not be relied upon.

In the event of a breach of the foregoing warranty, Millipore's sole obligation shall be to repair or replace, at its option, the applicable product or part thereof, provided the customer notifies Millipore promptly of any such breach. If after exercising reasonable efforts, Millipore is unable to repair or replace the product or part, then Millipore shall refund to the Company all monies paid for such applicable Product. **MILLIPORE SHALL NOT BE LIABLE FOR CONSEQUENTIAL, INCIDENTAL, SPECIAL OR ANY OTHER DAMAGES RESULTING FROM ECONOMIC LOSS OR PROPERTY DAMAGE SUSTAINED BY ANY COMPANY CUSTOMER FROM THE USE OF ITS PRODUCTS.**

(c) 2008: Millipore Corporation. All rights reserved. No part of these works may be reproduced in any form without permission in writing



