

SENSITIVE HUMAN LEPTIN RIA KIT
250 TUBES (Cat. # SHL-81K)

I.	Intended Use	2
II.	Principles Of Procedure	2
III.	Reagents Supplied	2
IV.	Storage and Stability	3
V.	Reagent Precautions	3
VI.	Materials Required But Not Provided	4
VII.	Specimen Collection And Storage	4
VIII.	Assay Procedure	5
IX.	Calculation	6
X.	Interpretation	6
XI.	Normal Fasting Range	6
XII.	Assay Characteristics	7
XIII.	Quality Controls	11
XIV.	Replacement Reagents	11
XV.	Ordering Information	12
XVI.	References	13
XVII.	Kits Available by Linco Research	13

**SENSITIVE HUMAN LEPTIN RIA KIT
250 TUBES (Cat. # SHL-81K)**

I. INTENDED USE

This Sensitive Human Leptin Radioimmunoassay (RIA) Kit utilizes an antibody made specifically against Human Leptin. Sensitivity of 0.05 ng/ml can be achieved when using a 100 μ l sample of serum, plasma, or tissue culture media, in a two-day, disequilibrium assay (400 μ l Total Volume). *This kit is for research purposes only.*

II. PRINCIPLES OF PROCEDURE

In radioimmunoassay, a fixed concentration of labeled tracer antigen is incubated with a constant dilution of antiserum such that the concentration of antigen binding sites on the antibody is limited, for example, only 50% of the total tracer concentration may be bound by antibody. If unlabeled antigen is added to this system, there is competition between labeled tracer and unlabeled antigen for the limited and constant number of binding sites on the antibody. Thus, the amount of tracer bound to antibody will decrease as the concentration of unlabeled antigen increases. This can be measured after separating antibody-bound from free tracer and counting one or the other, or both fractions. A calibration or standard curve is set up with increasing concentrations of standard unlabeled antigen and from this curve the amount of antigen in unknown samples can be calculated. Thus, the four basic necessities for a radioimmunoassay system are: a specific antiserum to the antigen to be measured, the availability of a radioactive labeled form of the antigen, a method whereby antibody-bound tracer can be separated from the unbound tracer, and finally, an instrument to count radioactivity.

The Linco Research Sensitive Human Leptin assay utilizes 125 I-labeled Human Leptin and a Sensitive Human Leptin antiserum to determine the level of Leptin in serum, plasma or tissue culture media by the double antibody/PEG technique.

III. REAGENTS SUPPLIED

Each kit is sufficient to run 250 tubes and contains the following reagents.

A. Assay Buffer

0.05M Phosphosaline pH 7.4 containing 0.025M EDTA, 0.08% Sodium Azide, 1% RIA Grade BSA and 0.05% Triton X-100
Quantity: 40 ml/vial
Preparation: Ready to use

B. Antiserum

Guinea Pig anti-Sensitive Human Leptin Serum in Assay Buffer
Quantity: 26 ml/vial
Preparation: Ready to use

C. 125 I-Human Leptin

125 I-Human Leptin Label, HPLC purified (specific activity 135 μ Ci/ μ g)
Lyophilized for stability. Freshly iodinated label contains <3 μ Ci, (<111 kBq) calibrated to the 1st Monday of each month.
Quantity: 27 ml/vial upon hydration
Preparation: Contents Lyophilized. Hydrate with entire contents of Label Hydrating Buffer. Allow to sit at room temperature for 30 minutes, with occasional gentle mixing.

D. Label Hydrating Buffer

Assay Buffer containing Normal Guinea Pig IgG as a carrier. Used to hydrate 125 I-Human Leptin
Quantity: 27 ml/vial
Preparation: Ready to use

III. REAGENTS SUPPLIED (continued)

E. Standards

Purified Recombinant Human Leptin in Assay Buffer at the following concentrations:

0.05, 0.1, 0.2, 0.5, 1.0, 2.0, 5.0, 10.0 ng/ml

Quantity: 1 ml/vial

Preparation: Ready to use

F. Quality Controls 1 & 2

Purified Recombinant Human Leptin in Assay Buffer

Quantity: 1 ml/vial

Preparation: Ready to use

G. Precipitating Reagent

Goat anti-Guinea Pig IgG serum, 3% PEG and 0.05% Triton X-100 in 0.05M Phosphosaline, 0.025M EDTA, 0.08% Sodium Azide

Quantity: 260 ml/vial

Preparation: Ready to use; chill to 4°C.

IV. STORAGE AND STABILITY

Refrigerate all reagents between 2 and 8°C for short-term storage. For prolonged storage (>2 weeks), freeze at $\leq -20^{\circ}\text{C}$. Avoid multiple (>5) freeze/thaw cycles. Refer to date on bottle for expiration when stored at $\leq -20^{\circ}\text{C}$. Do not mix reagents from different kits unless they have the same lot number.

V. REAGENT PRECAUTIONS

A. Radioactive Materials

This radioactive material may be received, acquired, possessed and used only by research personnel or clinical laboratories for in vitro research tests not involving internal or external administration of the material, or the radiation there from, to human beings or animals. Its receipt, acquisition, possession, use and transfer are subject to the regulations of the U. S. Nuclear Regulatory Commission (NRC) or of a State with which the Commission has entered into an agreement for the exercise of regulatory authority.

The following are suggested general rules for the safe use of radioactive material. The customer's Radiation Safety Officer (RSO) is ultimately responsible for the safe handling and use of radioactive material.

1. Wear appropriate personal devices at all times while in areas where radioactive materials are used or stored.
2. Wear laboratory coats, disposable gloves, and other protective clothing at all times.
3. Monitor hands, shoes, and clothing and immediate area surrounding the workstation for contamination after each procedure and before leaving the area.
4. Do not eat, drink, or smoke in any area where radioactive materials are stored or used.
5. Never pipette radioactive material by mouth.
6. Dispose of radioactive waste in accordance with NRC rules and regulations.
7. Avoid contaminating objects such as telephones, light switches, doorknobs, etc.
8. Use absorbent pads for containing and easily disposing of small amounts of contamination.
9. Wipe up all spills immediately and thoroughly and dispose of the contaminated materials as radioactive waste. Inform Radiation Safety Officer.

V. REAGENT PRECAUTIONS (continued)

B. Sodium Azide

Sodium Azide has been added to all reagents as a preservative at a concentration of 0.08%. Although it is at a minimum concentration, Sodium Azide may react with lead and copper plumbing to form highly explosive metal azides. On disposal, flush with a large volume of water to prevent azide build up.

VI. MATERIALS REQUIRED BUT NOT PROVIDED

1. Borosilicate glass tubes, 12 x 75 mm. (NOTE: Polypropylene or polystyrene tubes may be used if the investigator finds that the pellet formation is acceptably stable in their system.)
2. 100 µl pipet with disposable tips
3. 100 µl & 1.0 ml repeating dispenser
4. Refrigerated swing bucket centrifuge capable of developing 2,000 - 3,000 xg. (Use of fixed-angle buckets are not recommended.)
5. Absorbent paper
6. Vortex mixer
7. Refrigerator
8. Gamma Counter

VII. SPECIMEN COLLECTION AND STORAGE

1. A maximum of 100 µl per assay tube of serum or plasma can be used, although, 50 µl per assay tube is adequate for most applications. Tissue culture and other media may also be used.
2. Care must be taken when using heparin as an anticoagulant, since an excess will provide falsely high values. Use no more than 10 IU heparin per ml of blood collected.
3. Specimens can be stored at 4°C if they will be tested within 24 hours of collection. For longer storage, specimens should be stored at ≤ -20°C. Avoid multiple (>5) freeze/thaw cycles.
4. Avoid using samples with gross hemolysis or lipemia.

VIII. ASSAY PROCEDURE

For optimal results, accurate pipetting and adherence to the protocol are recommended.

Day One

1. Pipette 300 µl of Assay Buffer to the Non-Specific Binding (NSB) tubes (3-4). Pipette 200 µl of Assay Buffer in the Reference (Bo) tubes (5-6). Pipette 100 µl of Assay Buffer to tubes seven through the end of the assay.
2. Pipette 100 µl of Standards and Quality Controls in duplicate (see assay flow chart).
3. Pipette 100 µl of each sample in duplicate. (NOTE: Smaller volumes of sample may be used when Leptin concentrations are anticipated to be elevated or when sample size is limited. Additional Assay Buffer should be added to compensate for the difference so that the volume is equivalent to 100 µl (e.g., when using 50 µl of sample, add 50 µl of Assay Buffer). Refer to Section IX for calculation modification.
4. Pipette 100 µl of Sensitive Human Leptin Antibody to all tubes except Total Count tubes (1-2) and NSB tubes (3-4).
5. Vortex, cover, and incubate overnight (20-24 hours) at room temperature.

Day Two

6. Pipet 100 µl of ¹²⁵I-Human Leptin tracer to all tubes. Important: For preparation, see Section III, Part C.
7. Vortex, cover, and incubate overnight (22-24 hours) at room temperature.

Day Three

8. Add 1.0 ml of cold (4°C) Precipitating Reagent to all tubes except Total Count tubes (1-2).
9. Vortex and incubate 20 minutes at 4°C.
10. Centrifuge, at 4°C, for 20 minutes at 2,000-3,000 xg. Note: If less than 2,000 xg is used, the time of centrifugation must be increased to obtain a firm pellet (e.g. 40 minutes). Multiple centrifuge runs within an assay must be consistent. Conversion of rpm to xg:

$$xg = (1.12 \times 10^{-5}) (r) (rpm)^2$$
 r = radial distance in cm (from axis of rotation to the bottom of the tube)
 rpm = revolutions per minute
11. Immediately decant supernatant from all centrifuged tubes except Total Count tubes (1-2). Drain tubes for 15-60 seconds (be consistent between racks), and blot excess liquid from lip of tubes. Note: Invert tubes only one time. Pellets are fragile and slipping may occur.
12. Count all tubes in a gamma counter for 1 minute. Calculate the ng/ml of Sensitive Human Leptin in unknown samples using automated data reduction procedures (see Section IX).

Assay Procedure Flow Chart

Day One					Day Two		Day Three	
Set-up	Step 1	Step 2&3	Step 4	Step 5	Step 6	Step 7	Step 8	Steps 9-12
Tube Number	Add Assay Buffer	Add Standard/QC Sample	Add Sensitive Human Leptin Antibody	Vortex, Cover, and Incubate 20-24 hrs at RT	Add I-125 Human Leptin Tracer	Vortex, Cover and Incubate 22-24 hrs at RT	Add Precipitating Reagent	Incubate 20 min. at 4°C, Centrifuge at 4°C for 20 min Decant and Count
1,2	-	-	-		100 µl		-	
3,4	300 µl	-	-		100 µl		1.0 ml	
5,6	200 µl	-	100 µl		100 µl		1.0 ml	
7,8	100 µl	100 µl of 0.05 ng/ml	100 µl		100 µl		1.0 ml	
9,10	100 µl	100 µl of 0.1 ng/ml	100 µl		100 µl		1.0 ml	
11,12	100 µl	100 µl of 0.2 ng/ml	100 µl		100 µl		1.0 ml	
13,14	100 µl	100 µl of 0.5 ng/ml	100 µl		100 µl		1.0 ml	
15,16	100 µl	100 µl of 1.0 ng/ml	100 µl		100 µl		1.0 ml	
17,18	100 µl	100 µl of 2.0 ng/ml	100 µl		100 µl		1.0 ml	
19,20	100 µl	100 µl of 5.0 ng/ml	100 µl		100 µl		1.0 ml	
21,22	100 µl	100 µl of 10.0 ng/ml	100 µl		100 µl		1.0 ml	
23,24	100 µl	100 µl of QC 1	100 µl		100 µl		1.0 ml	
25,26	100 µl	100 µl of QC 2	100 µl		100 µl		1.0 ml	
27,28	100 µl	100 µl of unknown	100 µl		100 µl		1.0 ml	

IX. CALCULATIONS AND TRANSFORMATIONS

A. Explanation

The calculations for Sensitive Human Leptin can be automatically performed by most gamma counters possessing data reduction capabilities or by independent treatment of the raw data using a commercially available software package⁵. Choose weighted 4-parameter or weighted log/logit for the mathematical treatment of the data. [NOTE: Be certain the procedure used subtracts the NSB counts from each average count, except Total Counts, prior to final data reduction.]

B. Manual Calculation

1. Average duplicate counts for Total Count tubes (1-2), NSB tubes (3-4), Total Binding tubes (reference, Bo) (5-6), and all duplicate tubes for standards and samples to the end of the assay.
2. Subtract the average NSB counts from each average count (except for Total Counts). These counts are used in the following calculations.
3. Calculate the percentage of tracer bound
[(Total Binding Counts/Total Counts) X 100].
This should be 35-50%.
4. Calculate the percentage of total binding (%B/Bo) for each standard and sample
[%B/Bo = (Sample or Standard/Total Binding) X 100]
5. Plot the % B/Bo for each standard on the y-axis and the known concentration of the standard on the x-axis. The use of log-log graph paper will result in a nearly linear curve.
6. Construct the reference curve by joining the points with a smooth curve.
7. Determine the ng/ml of Sensitive Human Leptin in the unknown samples and controls by interpolation of the reference curve.

[NOTE: When sample volumes assayed differ from 100 µl, an appropriate mathematical adjustment must be made to accommodate for the dilution factor (e.g., if 50 µl of sample is used, then calculated data must be multiplied by 2).]

X. INTERPRETATION

A. Acceptance Criteria

1. The run will be considered accepted when all Quality Control values fall within the calculated Quality Control Range; if any QC's fall outside the control range, review results with the supervisor.
2. If the difference between duplicate results of a sample is >10% CV, repeat the sample.
3. The limit of sensitivity for the Sensitive Human Leptin assay is 0.05 ng/ml (100 µl sample size).
4. The limit of linearity for the Sensitive Human Leptin assay is 10.0 ng/ml (100 µl sample size). Any result greater than 10.0 ng/ml should be repeated on dilution using Assay Buffer as a diluent.

XI. NORMAL FASTING RANGE

Leptin levels are directly correlated with degree of adiposity.

Mean Leptin Values (BMI ranges 18-25):

Lean Men 3.8 ± 1.8 ng/ml

Lean Women 7.4 ± 3.7 ng/ml

Levels rise approximately 2.5 times faster in women per unit BMI as compared to men.

XII. ASSAY CHARACTERISTICS

A. Sensitivity

The lowest level of Leptin that can be detected by this assay is 0.05 ng/ml when using a 100 µl sample size.

B. Performance

The following parameters of assay performance are expressed as Mean ± Standard Deviation.

$$ED_{80} = 0.28 \pm 0.04 \text{ ng/ml}$$

$$ED_{50} = 1.05 \pm 0.12 \text{ ng/ml}$$

$$ED_{20} = 4.48 \pm 0.66 \text{ ng/ml}$$

C. Specificity

The specificity (also known as selectivity) of an analytical test is its ability to selectively measure the analyte in the presence of other like components in the sample matrix.

Human Leptin	100%
Rat Leptin	< 1%
Mouse Leptin	< 2%
Canine Leptin	< 0.2%
Human Insulin	*
Glucagon	*

*not detectable

D. Precision

Within and Between Assay Variation

Sample No.	Mean ng/ml	Within % CV	Between % CV
1	0.44	5.25	8.90
2	0.92	3.74	3.47
3	1.50	5.69	3.24
4	3.14	7.48	4.02
5	4.24	5.97	8.67

Within and between assay variation was performed on five human serum samples containing varying concentrations of Human Leptin. Data (mean and % CV) shown are from five duplicate determinations of each serum sample in five separate assays.

XII. ASSAY CHARACTERISTICS (continued)

E. Recovery

Spiking Recovery of Leptin in Human Serum

Sample No.	Leptin Added ng/ml	Observed ng/ml	Expected ng/ml	% Recovery
1	0	0.30	-	-
2	0.2	0.52	0.50	104
3	0.5	0.85	0.80	106
4	1.0	1.50	1.30	115
5	2.0	2.72	2.30	118

Varying concentrations of Human Leptin were added to five human serum samples and the leptin content was determined by RIA. Mean of the observed levels from five duplicate determinations in five separate assays are shown. Percent recovery was calculated on the observed vs. expected.

F. Linearity

Effect of Serum Dilution


Sample #	Volume Sampled	Observed ng/ml	Expected ng/ml	% Of Expectd
1	100 µl	1.07	1.07	100
	75 µl	0.99		93
	50 µl	0.88		82
	25 µl	0.78		73
2	100 µl	2.07	2.07	100
	75 µl	2.00		97
	50 µl	1.85		89
	25 µl	1.50		72
3	100 µl	3.14	3.14	100
	75 µl	3.06		97
	50 µl	3.05		97
	25 µl	2.54		81
4	100 µl	4.24	4.24	100
	75 µl	4.34		102
	50 µl	4.42		104
	25 µl	4.05		96

Aliquots of pooled human serum containing varying concentrations of Leptin were analyzed in the volumes indicated. Dilution factors of 1.0, 1.33, 2.0 and 4.0 representing 100 µl, 75 µl, 50 µl and 25 µl, respectively, were applied in calculating observed concentrations. Mean Leptin levels and percent of expected for five separate assays are shown.

XII. ASSAY CHARACTERISTICS (continued)

G. Correlation Between Linco Kits

Comparison of Human Serum Leptin Analyzed by Human Leptin RIA kit (Cat. # HL-81K) and Sensitive Human Leptin RIA kit (Cat. # SHL-81K).

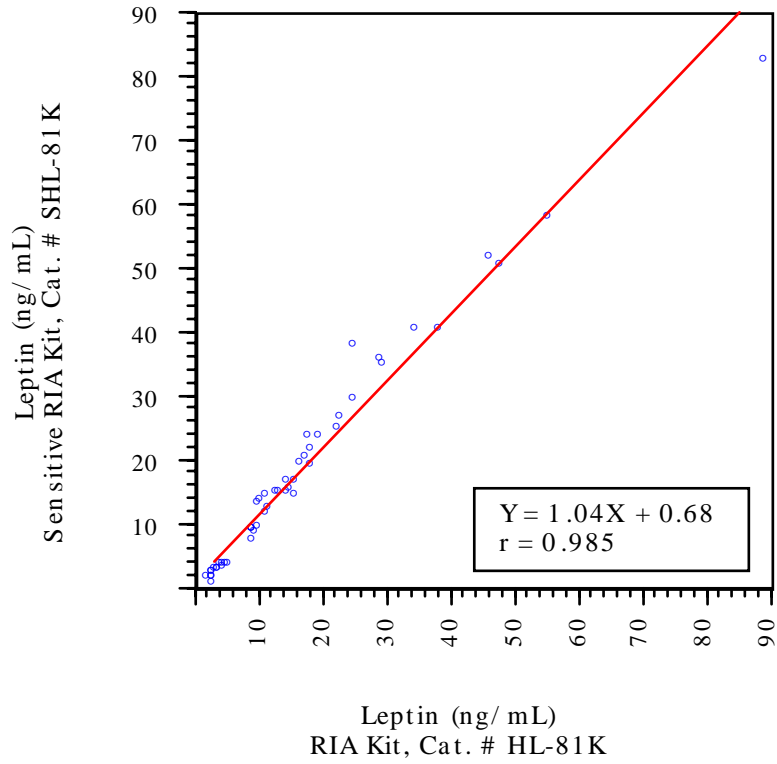

XII. ASSAY CHARACTERISTICS (continued)

H. Example of Assay Results

This data is presented as an example only and should not be used in lieu of a standard curve prepared with each assay.

Tube #	ID	CPM	Ave CPM	Ave Net CPM	% B/Bo	ng/ml
1	Totals	18253				
2	"	22772	20513			
3	NSB	618				
4	"	648	633			
5	Bo	9523				
6	"	9472	9498	8865		
<u>Standards</u>						
7	0.05 ng/ml	9134				
8		9011	9073	8440	95.2	
9	0.1 ng/ml	8777				
10		8704	8741	8108	91.5	
11	0.2 ng/ml	8220				
12		8227	8224	7591	85.6	
13	0.5 ng/ml	6692				
14		6664	6678	6045	68.2	
15	1.0 ng/ml	5024				
16		5135	5080	4447	50.2	
17	2.0 ng/ml	3390				
18		3310	3350	2717	30.7	
19	5.0 ng/ml	2325				
20		2236	2281	1648	18.6	
21	10.0 ng/ml	1908				
22		1904	1906	1273	14.4	
<u>Controls/Unknown</u>						
23	QC 1	7701				
24		7378	7540	6907	77.9	0.27
25	QC 2	3456				
26		3391	3424	2791	31.5	2.39
27-n	Unknown					

XIII. QUALITY CONTROLS

Good Laboratory Practice (GLP) requires that Quality Control (QC) specimens be run with each standard curve to check the assay performance. Two levels of controls are provided for this purpose. These and any other control materials should be assayed repeatedly to establish mean values and acceptable ranges. Each individual laboratory is responsible for defining their system for quality control decisions and is also responsible for making this system a written part of their laboratory manual.

Recommended batch analysis decision using two controls (Westgard Rules)⁶:

1. When both controls are within ± 2 SD. Decision: Approve batch and release analyte results.
2. When one control is outside ± 2 SD and the second control is within ± 2 SD.
Decision: Hold results, check with supervisor. If no obvious source of error is identified by the below mentioned check of systems, the supervisor may decide to release the results.

Technician check of systems:

1. Check for calculation errors
2. Repeat standards and controls
3. Check reagent solutions
4. Check instrument

XIV. REPLACEMENT REAGENTS

Reagents	Cat #
¹²⁵ I-Human Leptin (<3 μ Ci, <111 kBq)	9081
Label Hydrating Buffer (27 ml)	LHB-P
Human Leptin Standards (1 ml each)	8081S-K
GP anti-Human Leptin (Sensitive) Antibody (26 ml)	1081S-K
Precipitating Reagent (260 ml)	PR-UV
Quality Control 1 & 2 (1 ml each)	6000S-K
Assay Buffer (40 ml)	AB-PTR

XV. ORDERING INFORMATION

A. To place an order:

For USA Customers:

Please provide the following information to our customer service department to expedite your telephone, fax or mail order:

1. Your name, telephone and/or fax number
2. Customer account number
3. Shipping and billing address
4. Purchase order number
5. Catalog number and description of product
6. Quantity and product size

NOTE: Appropriate license from NRC (or equivalent) must be on file at LINCO before radioactive orders can be shipped.

TELEPHONE ORDERS:

Toll free US (866) 441-8400
(636) 441-8400

FAX ORDERS: (636) 441-8050

MAIL ORDERS: LINCO Research
6 Research Park Drive
St. Charles, Missouri 63304 U.S.A.

For International Customers:

To best serve our international customers, it is LINCO's policy to sell our products only through a network of distributors. To place an order or to obtain additional information about LINCO products, please contact your local distributor.

B. Conditions of Sale

All products are for research or manufacturing use only. They are not intended for use in clinical diagnosis or for administration to humans or animals. All products are intended for *in vitro* use only.

C. Terms

Invoices are net 30 days in U.S. dollars. Prices are subject to change without notice.

D. Material Safety Data Sheets (MSDS)

Material safety data sheets for LINCO Research products may be ordered by fax or phone. See section A above for details on ordering.

XVI. REFERENCES

1. Pelleymounter, M.A., et. al. Effects of the Obese Gene Product on Body Weight Regulation in Ob/ob Mice. *Science*. 269:540-543, 1995.
2. Maffei, M., et. al. Leptin Levels in Human and Rodent: Measurement of Plasma Leptin and ob RNA in Obese and Weight-reduced Subjects. *Nature Med.* Vol. 1, 11:1155-1611, 1995.
3. Ma, Zhongmin, et al. Radioimmunoassay of Leptin in Human Plasma. *Clinical Chemistry*. 42:942-946, June, 1996.
4. Thorell, J.I. *Scand. J. Clin. Lab. Invest.* 31:187, 1973.
5. Feldman, H. and Rodbard, D. "Mathematical Theory of Radioimmunoassay", in: W.D Odell and Doughaday, W.H. (Ed.), Principles of Competitive Protein-Binding Assays. Philadelphia: J.B. Leppincott Company; pp 158-203, 1971.
6. Westgard, J.O., et. al. A Multi-rule Shewhart Chart for Quality Control in Clinical Chemistry. *Clin. Chem.* 27:493-501, 1981.

XVII. KITS AVAILABLE FROM LINCO RESEARCH

Radioimmunoassays:

Human Leptin	Ultra Sensitive Human Insulin
Sensitive Human Leptin	Human Proinsulin
Mouse Leptin	Canine C-Peptide
Multi Species Leptin	Human C-Peptide
Rat Leptin	Porcine C-Peptide
Primate Leptin	Rat C-Peptide
Rat Insulin	Glucagon
Sensitive Rat Insulin	Glucagon Like Peptide-1 (Total)
Porcine Insulin	Glucagon Like Peptide-1 (Active)
Lispro Insulin	Rat Growth Hormone
Human Insulin Specific	Porcine/Canine Growth Hormone
Mouse Adiponectin	Human Adiponectin
Human Ghrelin (Total)	Human Ghrelin (Active)
Mouse Resistin	Human PYY (Total)

ELISAs:

Human Amylin (Active)	Human Leptin
Human Amylin (Total)	Glucagon Like Peptide-1 (Active)
Human Insulin	Human C-Peptide
Rat/Mouse Insulin	Human Resistin
Mouse Adiponectin	Human Adiponectin
Human Proinsulin (Total)	Human Proinsulin (Intact)
Desacyl-Ghrelin	Active Ghrelin

XVII. KITS AVAILABLE FROM LINCO RESEARCH (continued)

Multiplex:

Human Cytokine (IL-1 α , IL-1 β , IL-2, IL-4, IL-5, IL-6, IL-7, IL-8, IL-10, IL-12(p70), IL-13, IL-15, IL-17, IFN γ , GM-CSF, TNF α , IP-10, G-CSF, Eotaxin, MCP-1, MIP-1 α , RANTES), Premixed

Mouse Cytokine (IL-1 α , IL-1 β , IL-2, IL-4, IL-5, IL-6, IL-7, IL-9, IL-10, IL-12(p70), IL-13, IL-15, IL-17, IFN γ , IP-10, G-CSF, GMCSF, TNF α , MIP-1 α , MCP-1, KC, RANTES), Premixed

Rat Cytokine (IL-1 β , IL-4, IL-6, IL-10, IFN γ , GRO-KC, TNF α , GM-CSF, IL-1 α , MCP-1, IL-2, IL-12, IL-5, IL-18), Premixed

Human Endocrine (Leptin, GLP-1, C-Peptide, Insulin, Glucagon, Amylin (Total), Amylin (Active))

Rat Endocrine (Insulin, Amylin, Leptin, Glucagon, GLP-1,)

Mouse Endocrine (Insulin, Amylin, Leptin, Glucagon, GLP-1)

Apolipoproteins (Apo AI, Apo AII, Apo B, Apo CII, Apo CIII, Apo E)

Mouse Adipokine (Insulin, MCP-1, Leptin, IL-6, TNF α , tPAI-1, Resistin)

Mouse Adipocyte (MCP-1, Leptin, IL-6, TNF α , tPAI-1, Resistin, Adiponectin)

Single plex:

Human Cytokine (IL-10, TNF α , RANTES)

Mouse Cytokine (TNF α)

Mouse Adipokine (Adiponectin)